

## notes and applications from Bioanalytical Systems, Inc.

2701 Kent Ave • W. Lafayette, IN 47906 • (765) 463-4527 • FAX (765) 497-1102 • [bas@bioanalytical.com](mailto:bas@bioanalytical.com)

The NDA should always be added last to minimize side product formation.

The reaction mixture should be protected from light as it can undergo photodegradation.

#### References

1. P. de Montigny, J.F. Stobaugh, R.S. Givens, R.G. Carlson, K. Srinivasachar, L.A. Stemson and T. Higuchi, *Anal. Chem.* 59 (1987) 1096-1101.
2. B.K. Matuszewski, R.S. Givens, K. Srinivasachar, R.G. Carlson and T. Higuchi, *Anal. Chem.* 59 (1987) 1102-1105.
3. S.M. Lunte and O.S. Wong, *Current Separations* 10 (1990) 19-26.

\* US PATENT #4837166

